

HISTOSTAR EMBEDDING STATION

Quick start guide

Please keep the following default settings:

“STATUS”

- **timer** always OFF
- **paraffin** always ON
- **hot spot** OFF when not in use
- **tissue tank** always ON
- **mold storage** OFF when not in use
- cooling option-> **cold plate** OFF when not in use
- cooling option-> **cold spot** OFF when not in use

5-10 min before you start

- switch ON cold plate and cold spot (cooling options)
- switch ON hot spot and mold storage

when you are done

- switch OFF cold plate and cold spot (cooling options)
- switch OFF hot spot and mold storage
- clean holds from paraffin and put into mold storage
- empty paraffin spill tray

Please do not switch off the instrument!